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12 Ihre Notizen

table of contents

1	introduction.....	26
2	certification of <i>piezosystem jena</i>	26
3	declaration of conformity.....	26
4	purchased part package.....	27
5	instructions for using piezo electrical elements and power supplies.....	27
6	safety instructions.....	28
6.1	installation, power supply.....	29
6.2	operation.....	29
6.3	maintenance and inspection.....	30
6.4	environmental conditions.....	30
7	instructions for checking the function of the system / quick start.....	30
8	how to operate the digital amplifier <i>NV100/D_NET</i>	31
8.1	common introduction.....	31
8.2	connections.....	31
8.3	function.....	32
8.4	technical data.....	35
8.5	pinning.....	36
8.6	communication.....	37
8.6.1	USB-communication.....	37
8.6.2	Ethernet-communication.....	37
8.6.3	Ethernet-configuration.....	38
8.7	commands.....	41
8.8	status register.....	42
8.9	error messages.....	42
9	handling.....	43
10	controller adjustment.....	44
11	troubleshooting.....	45
12	your notes.....	46

1 introduction

This manual describes the digital piezo amplifier series **NV100/D_NET** from **piezosystem jena**. You will also find additional information regarding piezoelectric products.

Definition: All systems from **piezosystem jena** such as electronics, actuators, and optical systems are called “units”.

If you have any problems please contact the manufacturer of the system: **piezosystem jena**, Stockholmer Str. 12, 07747 Jena. Phone: +49 36 41 66 88-0

2 certification of **piezosystem jena**

The company **piezosystem jena GmbH** has worked according to a DIN EN ISO 9001 certified quality management system since 1999. Its effectiveness is verified and proven by periodic audits by the TÜV.

This instruction manual includes important information for using piezo actuators and electronics. Please take the time to read this information. Piezo positioning systems are mechanical systems that offer the highest precision. Correct handling guarantees that this precision will be maintained over a long period of time.

3 declaration of conformity

The CE Declaration of Conformity is available on our website **www.piezosystem.com**.

4 purchased part package

- Please check the completeness of the delivery after receiving the shipment:
- piezo amplifier **NV100/D_NET**
- wide range power supply 24 V_{DC}
- USB cable
- instruction manual

For optional needed drivers or available Software see our homepage

5 instructions for using piezo electrical elements and power supplies

- Piezoelectric actuators from **piezosystem jena** are controlled by voltages up to 130V. These values can be quite hazardous. Therefore, read the installation instructions carefully and ensure that only authorized personnel handle the power supply.
- After transportation, piezoelectric actuators should be allowed to adapt to room temperature for approximately 2 hours before being switched on.
- Piezoelectric actuators are made from ceramic materials with and without metallic casings. The piezo-ceramic is a relatively brittle material. This should be noted when handling piezoelectric actuators. All piezo elements are sensitive to bending or shock forces.
- Due to the piezoelectric effect, piezo actuators can generate electrical charges by changing the mechanical load or the temperature, or by actions such as the ones described above.
- Piezoelectric actuators are able to work under high compressive forces. Only actuators with a pre-load can be used under tensile loads (these tensile forces must be less than the pre-load given in the data sheet). Please note that acceleration of the ceramic material (e.g., caused by fall down, discharging, or high dynamic application) will occur.
- After excitation of the actuators by a voltage in the upper control range, the ceramic will move and generate an opposite high voltage after disconnection.
- Heating of the ceramic material will occur during dynamic operation and is caused by structure conditional loss processes. This may cause failure if the temperature exceeds specified values cited below. With increasing temperature up to the Curie temperature T_C , (usual values approx. 140 °C – 250 °C) the piezoelectric effect disappears. We recommend working in temperatures up to $T_C/2$ (normally up to 80 °C).
- Piezoelectric actuators, such as stacks or other devices, work electrically as a capacitor. These elements are able to store electrical energy over a long period of time (up to some days) and the stored energy may be dangerous.
- If the actuator remains connected to the drive electronics, it will be unloaded within a second after shutdown and quickly reaches harmless voltage values.
- Piezo actuators can only generate voltages by warming or cooling (caused by the longitudinal change). The discharge potential should not be ignored due to the inner capacitance. This effect is insignificant at usual room temperature.
- Piezo actuators from **piezosystem jena** are adjusted and glued. Any opening of the unit will cause misalignment or possible malfunction and will result in the loss of the guarantee.
- Please only use original parts from **piezosystem jena**.

Please contact **piezosystem jena** or your local representative if there are any problems with your actuator or power supply.

Caution! Shock forces may damage the built-in ceramic elements. Please avoid such forces, and handle the units with care, otherwise the guarantee will be lost.

6 safety instructions

Icons:

RISK OF ELECTRIC SHOCK! Indicates that a risk of electric shock is present and the associated warning should be observed.

CAUTION! REFER TO OPERATOR'S MANUAL – Refer to your operator's manual for additional information, such as important operating and maintenance instructions.

RISK OF ELECTRIC SHOCK!

- Do not open the units! There are no user serviceable parts inside and opening or removing covers may expose you to dangerous shock hazards or other risks. Refer all servicing to qualified service personnel.
- Do not spill any liquids into the cabinet or use the units near water.

CAUTION!

- Allow adequate ventilation around the units so that heat can properly dissipate. Do not block ventilated openings or place the units near a radiator, oven, or other heat sources. Do not put anything on top of the units except those that are designed for that purpose (e.g. actuators).
- Only work with the units in a clean and dry environment! Only specially prepared units (e.g. actuators) can work under other conditions!
- Please only use original parts from **piezosystem jena**. **piezosystem jena** does not give any warranty for damages or malfunction caused by additional parts not supplied by **piezosystem jena**. Additional cables or connectors will change the calibration and other specified data. This can change the specified properties of the units and cause them to malfunction.
- Piezo elements are sensitive systems capable of the highest positioning accuracy. They will demonstrate their excellent properties only if they are handled correctly! Please mount them properly at the special mounting points.

Immediately unplug your unit from the wall outlet and refer servicing to qualified service personnel under the following conditions:

- when the cords or plugs are damaged
- if liquid has been spilled or objects have fallen into the unit
- if the unit has been exposed to rain or water
- if the unit has been dropped or the housing is damaged

6.1 installation, power supply

RISK OF ELECTRIC SHOCK

- Do not insert or unplug the power plug with wet hands, as this may result in electrical shock.
- Do not install in rooms where inflammable substances are stored. If flammable substances come into contact with electrical parts inside, it could result in fire or electrical shock.
- Do not damage or modify the power cord. Also, do not place heavy objects on the power cord, or pull on or excessively bend it, as this could cause electrical damage and result in a fire or electrical shock.
- Always grasp the plug portion when unplugging the power cord. Pulling on the power cord may expose or snap the core wire, or otherwise damage the power cord. If the cord is damaged, this could cause an electricity leak and result in a fire or electrical shock.

CAUTION!

- Do not use accessories other than the ones provided (e.g. power cord). Only plug the power cord into grounded power equipment and sockets.
- Do not place heavy objects on any cables (e.g. power cords, sensor cables, actuator cables, optical cables).
- Do not block ventilated openings or place the units near a radiator, oven, or other heat sources.
- Plug in the power cord completely so that it cannot loosen inadvertently.
- Leave sufficient space around the power plug so that it can be unplugged easily. If objects are placed around the power plug, you will be unable to unplug it in an emergency.
- Install the system so that the on/off-switch is easily accessible at all times.
- The power plug is the cut-off point to the main power supply.

6.2 operation

RISK OF ELECTRIC SHOCK!

- Do not open the units! There are no user serviceable parts inside and opening or removing covers may expose you to dangerous shock hazards or other risks. Refer all servicing to qualified service personnel.
- Do not spill inflammable substances inside the voltage amplifier. If these items come into contact with an electrical component inside the voltage amplifier, this may result in a fire or electrical shock.

CAUTION!

- If the voltage amplifier emits smoke, high heat, or unusual smells, immediately turn off the power switch and unplug the power plug from the outlet. Then contact our technical service.

6.3 maintenance and inspection

CAUTION! 

- Before cleaning the exterior box of the voltage amplifier, turn off the power switch and unplug the power plug. Failure to do so may result in a fire or electrical shock.
- Clean the exterior box using a damp cloth that has been firmly wrung-out. Do not use alcohols, benzene, paint thinner or other inflammable substances. If flammable substances come into contact with an electrical component inside the voltage amplifier, this may result in a fire or electrical shock.

6.4 environmental conditions

The amplifier can be used:

- indoors only
- at an altitude of up to 2000 m
- at a temperature between: 5 ... 35 °C
- at a relative humidity between: 5 ... 95 % (non-condensing)

The recommended environmental conditions:

- indoors only
- at an altitude of up to 2000 m
- at a temperature between: 20 ... 22 °C
- at a relative humidity between: 5 ... 80 % (non-condensing)

7 instructions for checking the function of the system / quick start

When you open the package, please check to make sure all the necessary parts are included (see packing list) and nothing is damaged. Check the electronics and the actuator for any visible damage:

- The top and bottom plates of the actuator (if it does not have another shape) should be parallel each to each other and not have any scratches.
- If there is any damage to the system please contact our local representative immediately!
- If the packaging material is damaged please confirm this with the shipping company.
- Connect the piezo actuator by using the D-SUB 15pin connector. Be sure the cables are connected properly to the electronics.
- Connect the device to PC with an USB cable.
- Connect the power supply (24 V_{DC})
- Normally, the PC installs necessary drivers automatically
After initialization, the LED lights green (open loop mode) or orange (closed loop mode)
- In case of irregularities, please read the chapter on *troubleshooting*.
- The function check is complete.

8 how to operate the digital amplifier *NV100/D_NET*


8.1 common introduction

The digital amplifier *NV100/D_NET* has designed to control the D-actuators from **piezosystem jena**. it can be used with actuators with or without integrated position sensor. Actuators with integrated position sensor allows positioning without creeping or hysteresis by digital PID loop control. The amplifier is controlled by USB only and reach a resolution of 16-bit. piezosystem jena supports driver for Windows® operating systems.

The integrated digital loop controller comes as PID controller, additionally a slew rate limiter and a low pass filter 4th order for set point filtering are available.


The digital amplifier *NV100/D_NET* allows the interchangeability from D-actuators and amplifier. All needed information is stored on the ID-chip, located in actuator connector, and will read by piezo controller after system start.

8.2 connections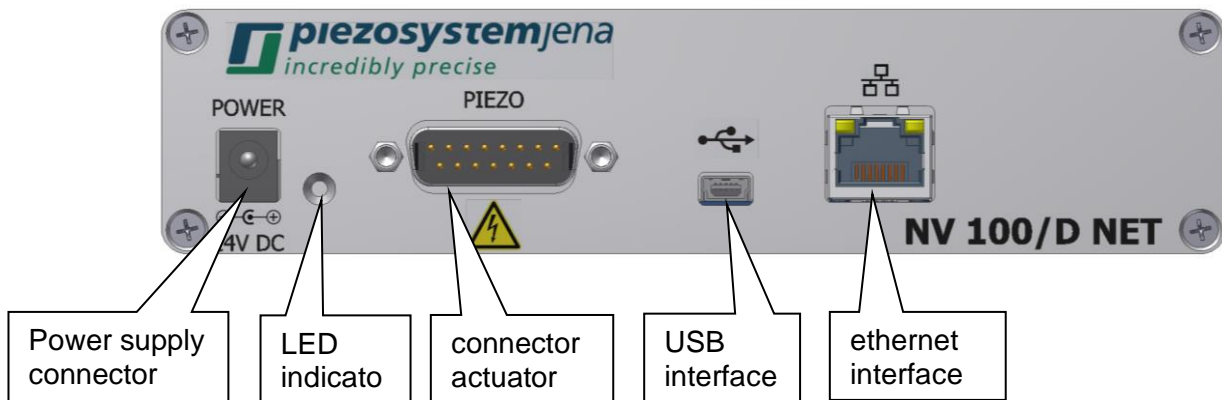

8.3 function

The following sketch shows the function groups of the amplifier: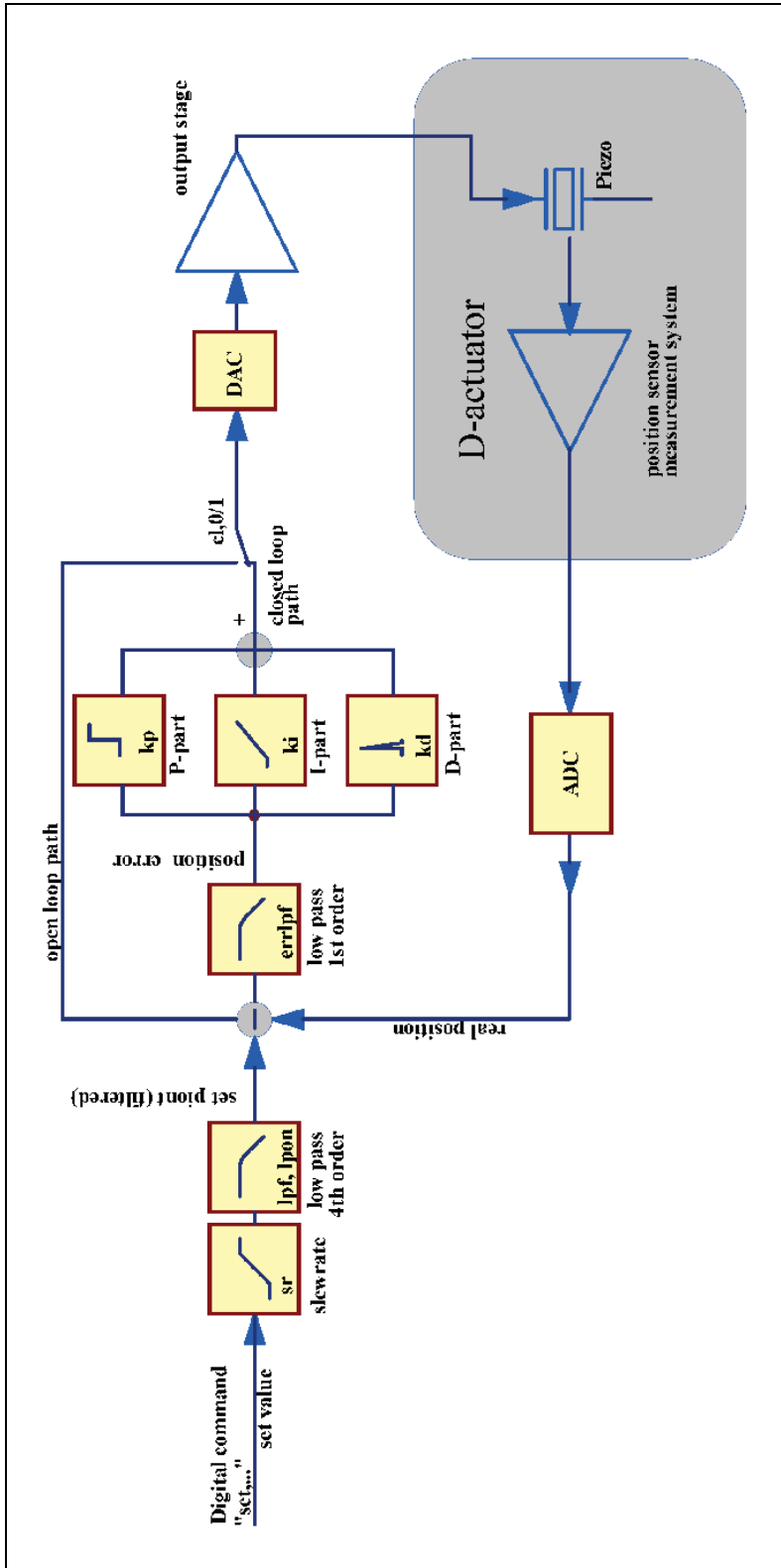

sketch 1: function groups of the amplifier

The controller can work in two different modes: The open loop mode (OL) and the closed loop mode (CL). In open loop mode, the piezo actuator is controlled by a voltage between -20 and +130 Volt. The filters “slew rate” and “low pass” can be used. The resulting stroke depends on the characteristic of piezo, it contains creeping and hysteresis.

In closed loop mode, these effects will be compensated by the digital loop controller. The functional groups for cl mode will be explained now:

The setpoint is given by USB command. The raise time can be adapted by slew rate limit and / or using the low pass filter 4th order to the mechanical System. These filters can be used to avoid excitation of oscillations before the signal reaches the control loop.

Now to the PID loop controller:

The PID-controller calculates the difference (err) between this input (com) and the value from the measurement system (mes).

$$\text{err} = \text{com} - \text{meas}$$

com = command
meas = position signal

The proportional term (kp) amplifies the control deviation without reducing frequency and drives the controller until the difference between the sensed motion and the command:

$$\text{yp} = \text{kp} * \text{err}$$

kp = adjustable gain
yp = output of p-term

The P-controller can't eliminate the error completely, because it needs a deviation to drive. To minimize the permanent offset of the P-term an integral action (I-term) is required. Ts is the time constant governing the time it takes for the output to get a certain value. For a step input it is the time taken for the output to equal the input.

$$\text{yi} = \text{ki} * \text{err} * \text{Ts}^2$$

Ts = 1 / sample frequency (20 kHz)
1/ki = time constant
yi = output of i-term
2= correction factor for compatibility to older devices

In general, the differential term combats oscillation (it adds damping) and increases the reaction speed of the PI-controller:

$$\text{yd} = \text{kd} * 1/\text{Ts} * (\text{err} - \text{err}[\text{n}-1])$$

err[n-1] = control deviation previous sample
kd = differential time constant
yd = output of d-term

Be careful when using the d-term, because stochastic errors, like noise, cause extreme reactions. The addition of these 3 terms generates a PID-controller. The output of the controller is the set value for the power amplifier to drive the piezo actuator:

$$\text{y} = \text{yp} + \text{yi} + \text{yd}$$

y = output of PID-controller


In open loop mode, the PID-controller is bridged and the command input controls the amplifier.

Using a subsequent steep notch filter the main resonant frequency will be damped by another order of magnitude. While using open loop you can also switch on the notch filter to reduce the stimulation of the resonant frequency.

The PID-specific parameters k_p , k_i , and k_d are stored in the actuator's ID-chip by default. These parameters will work for most applications. Depending on your special application and load, the mechanical response can be optimized.

The following sketch describes the step response of the different terms:

jump response of loop-controllers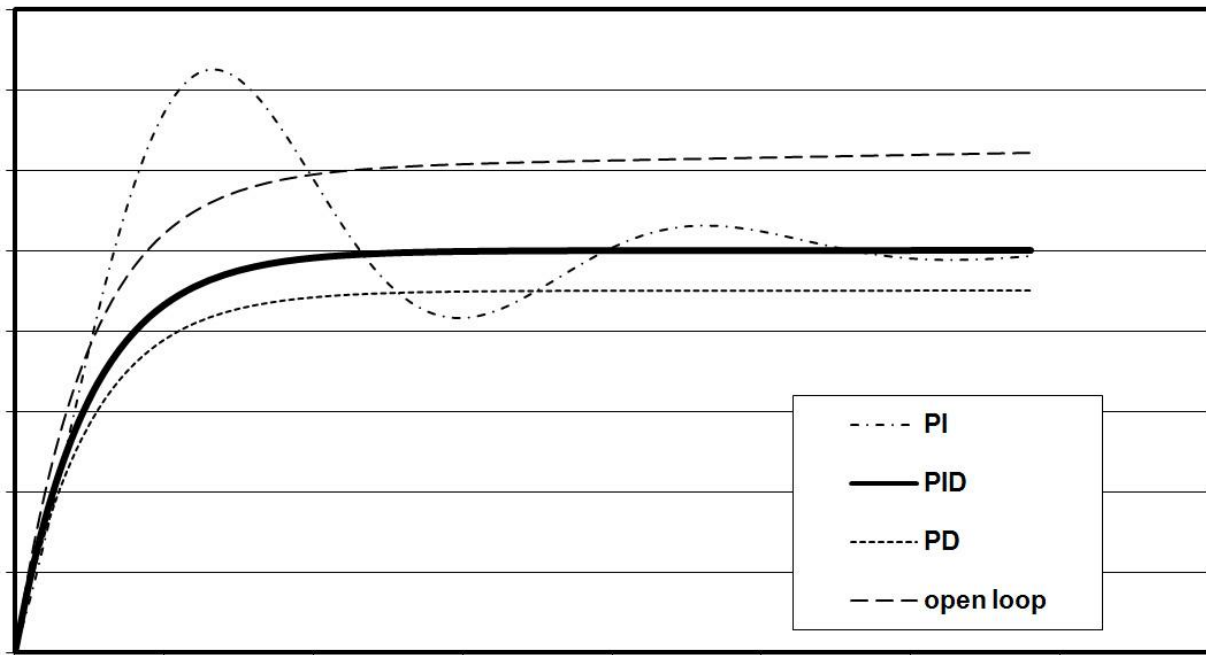

sketch 2: transfer function

The PID-specific parameters k_p , k_i , k_d and the several filters are stored in the actuator's ID-chip by default. These parameters will work for most applications. Depending on your special application and load, the mechanical response can be optimized.

8.4 technical data

	<i>NV100/D_NET</i>
Input voltage:	24 V _{DC}
Power supply connector:	low voltage socket with 2.1 mm pin
Input current:	max. 0.3 A @ 24 V, idle current 90 mA
Power consumption:	max. 8 W
Output power:	6.5 W (max. 12 W for NanoX® actuators)
Output voltages:	-20 V ... +130 V
Output current (constant current):	100 mA maximum (2x 50 mA for NanoX® actuators)
Output noise:	0.7 mV _{RMS} @ 500 Hz bandwidth
Actuator / measurement connector „PIEZO“:	15pin D-Sub plug
Safety functions:	short circuit proof, temperature fuse
Display (LED):	green / orange / red
Interface	USB, Ethernet
Interface connector	mini USB, RJ45
Cooling	passive
Dimensions (W H D):	165 x 38 x 120 mm ³
Weight:	0.35 kg

table: technical data

8.5 pinning

PIEZO: plug 15pin SUB-D

pin	designation	description
1, 2, 11	AGND	analog ground
3	+15V*	operating voltage measurement +15 V
4, 14	GND	digital ground
5	SDA	I ² C-Bus SDA
6	3.3V _{EEPROM} *	operating voltage ID-Chip
7	Vout2	for NanoX [®] actuator voltage 2 +130 ... -20 V
8	Vout	actuator voltage -20 ... +130 V
9	+MEAS	position signal -8 ... +8 V
10	-15V*	operating voltage measurement -15 V
12	SCL	I ² C-Bus SCL
13	DETECT	actuator detection
15	Piezo ground	actuator ground

[*] Not for external use


table: pinning

The actuator is connected to this plug. The actuator voltage goes to the piezo, the position signal from the measurement system goes to the DSP, and the integrated actuator information is stored on an ID-chip which goes to the main controller. It is important that the locking screws are screwed into the front panel. The voltages will be switched after connecting the actuator. When disconnecting the actuator, the dangerous Voltages will be switched off by a safety relay.

8.6 communication

8.6.1 USB-communication


The **NV100/D_NET** can easily communicate with your PC with the use of a terminal program such as HTerm. A USB connection cable (included) is required to connect to your computer. The properties of the COM port are: 115200 baud, 8 bit, no parity, 1 stop bit, software handshake (XON / XOFF).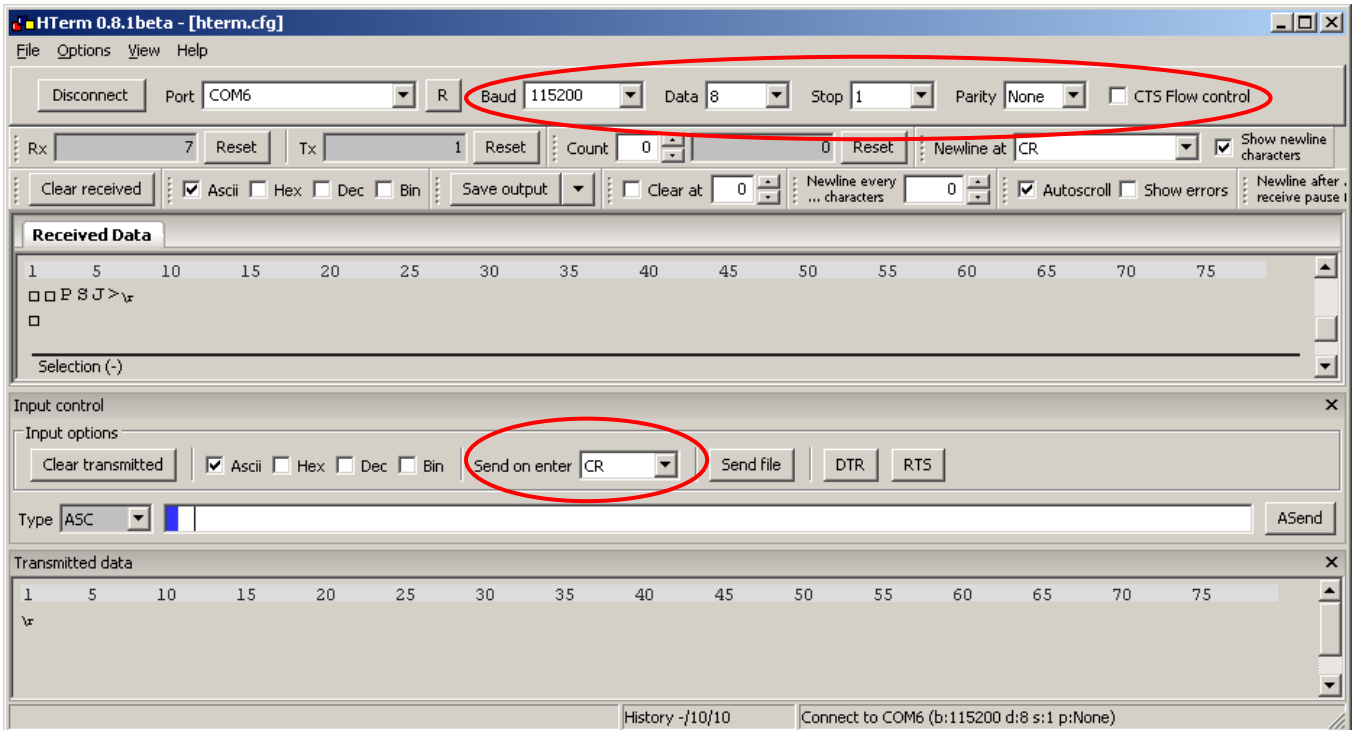


When the device is powered and connected, it answers after receiving “CR - carriage return” (Enter Key) with its prompt “NV100/D_NET>”

8.6.2 Ethernet-communication

The communication with the **NV100/D_NET** over ethernet can done with Windows own telnet. Alternative you can use other programmes like Putty. After the connection all commandos are possible like over USB. Default port is 23.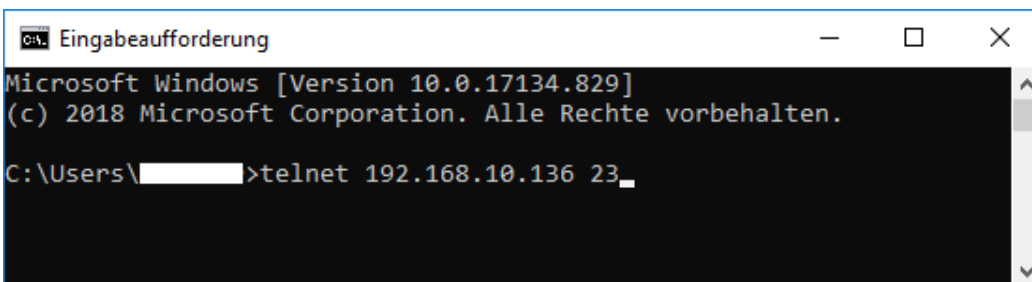


example: Telnet

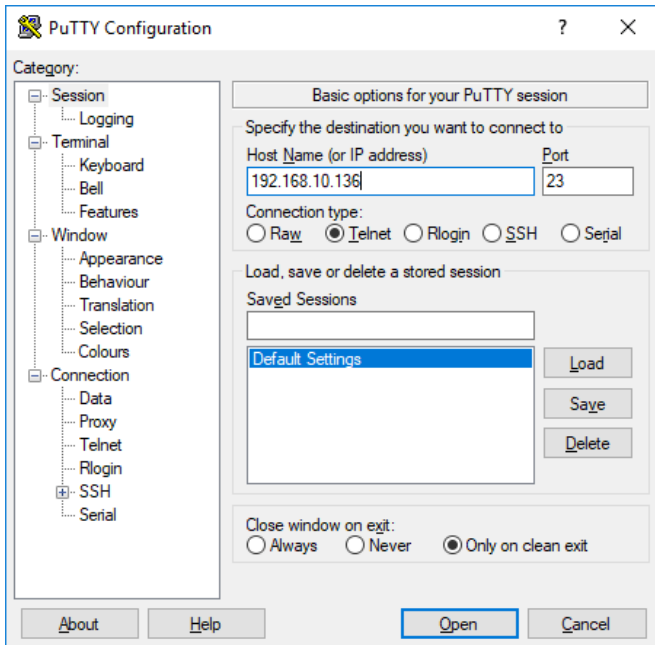example: Putty (Please change the ip-address)

The ip-address is dynamic over DHCP or can set static over the configuration. To get the ip-address of the device ask you network administrator or use our tool „IP-Search“. At problems with the firewall ask your network administrator.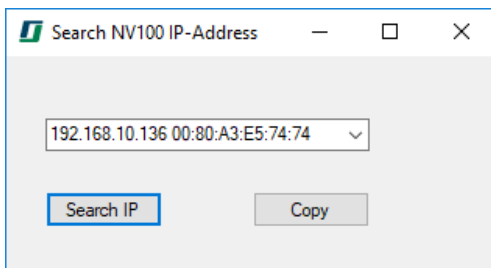


After the search of the IP-address you can copy it to clipboard with the „Copy“-button. They can later use in the command line for telnet.

Note: It is possible to use the host name, if the „DHCP FQDN option“ is enabled. If you don't set a name, the default is „C“ + the last 6 digit of the MAC address. (Example: MAC=00:80:A3:E5:74:74 → host name: „CE57474“)

8.6.3 Ethernet-configuration

The network setting can change over telnet (port: 9999) or with browser. There you can change the IP-address and the default port(23).

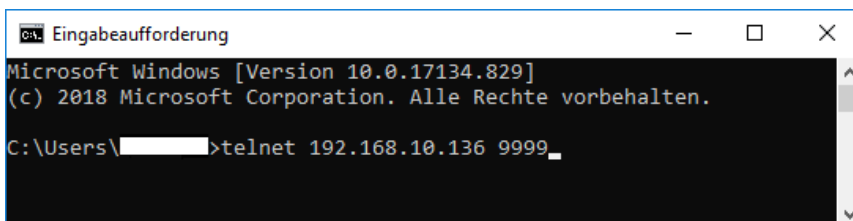In the first line you can read the MAC-address of the device. On the end you see the configuration menu.


```
Telnet 192.168.10.136
MAC address 0080A3E57474
Software version V6.10.0.3 (171229) XPTEXE

Press Enter for Setup Mode

*** basic parameters
Hardware: Ethernet TPI
IP addr - 0.0.0.0/DHCP/BOOTP/AutoIP, no gateway set
DNS Server not set
DHCP device name : CE57474
DHCP FQDN option: Disabled

*** Security
SNMP is          enabled
SNMP Community Name: public
Telnet Setup is  enabled
TFTP Download is enabled
Port 77FEh is    enabled
77FEh Access Mode is Read & Write
Web Server is    enabled
Web Setup is     enabled
ECHO is          disabled
Enhanced Password is disabled
Port 77F0h is    enabled

*** Channel 1
Baudrate 115200, I/F Mode 4C, Flow 00
Port 00023
Connect Mode : C0
Send '+++' in Modem Mode enabled
Show IP addr after 'RING' enabled
Auto increment source port disabled
Remote IP Adr: --- none ---, Port 00000
Disconn Mode : 00
Flush Mode : 00

*** Expert
TCP Keepalive   : 45s
ARP cache timeout: 600s
CPU performance: Regular
Monitor Mode @ bootup : enabled
RS485 tx enable : active low
HTTP Port Number : 80
SMTP Port Number : 25
MTU Size: 1400
TCP Re-transmission timeout: 500 ms
Alternate MAC: disabled
Ethernet connection type: auto-negotiate

*** E-mail
Mail server: 0.0.0.0
Unit       :
Domain     :
Recipient 1:
Recipient 2:

- Trigger 1
Serial trigger input: disabled
Channel: 1
Match: 00,00
Trigger input1: X
```

```
Telnet 192.168.10.136
Change Setup:
0 Server
1 Channel 1
3 E-mail
5 Expert
6 Security
7 Defaults
8 Exit without save
9 Save and exit          Your choice ?
```

With „Server“ you can set the IP-address to static. With „Channel 1“ you can change the default port (23).

With the ip address you can change the network settings in every browser.

CAUTION: In delivery condition the IP address is get over DHCP. In nets without DHCP server there is no access possible without change to a static ip. Please ask your network administrator.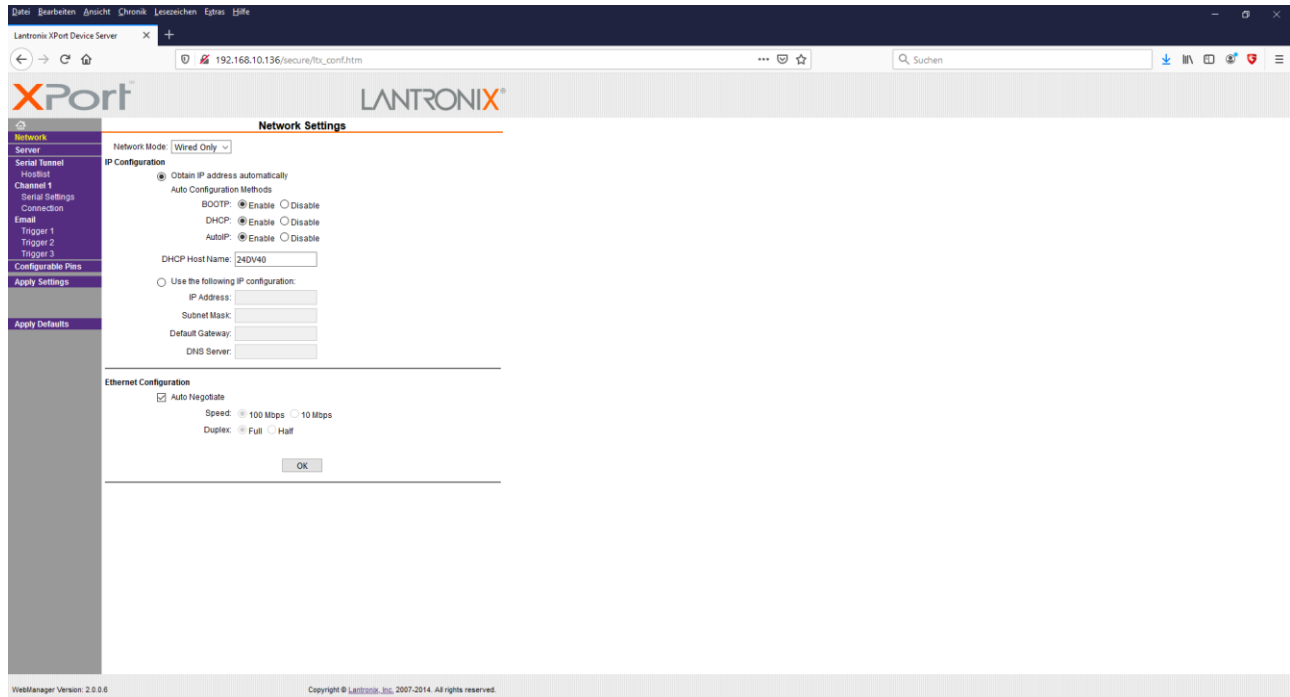

8.7 commands

commands with values: <command>,<value> Enter

command	description	value range	auto set after switch on
fenable	reformatting the actuator (soft start) after switch on	0= soft start disabled 1= soft start enabled	x
sinit	initial actuator position after switch on	0 ... 100%	x
set	command value: actuator voltage (ol) displacement (cl)	-20 to +130.000 [V] 0 to xxx.xxx [µm] (maximum actuator displacement, see datasheet)	-
cl	open loop / closed loop mode	0 = open loop 1 = closed loop	x
sr	slew rate limit	0.0000008...2000.0 [%/ms]	x
kp	proportional term	0 ... 10000	x
ki	integral term	0 ... 10000	x
kd	differential term	0 ... 10000	x
lpon	low pass filter for setpoint activating	0 = off 1 = on	x
lpf	low pass cut frequency	1 ... 10000 [Hz]	x
meas	position value	V (in ol mode) / actuator unit (in CL mode)	-
stat	request of device state	-	-
s	request of commands list	-	-

table: commands

8.8 status register

The status register is a 16bit register, in which each bit describes different properties of the amplifier or actuator. The decimal sum of all bits is the value of the status register, it can be requested by the “stat”- command.

bit	value	description	value decimal
0	2^0	0 – actuator not connected 1 - actuator connected	0 1
2,1	$2^2, 2^1$	0, 0 – actuator without position sensor 0, 1 – strain gauge sensor 1, 0 – capacitive sensor	0 2 4
3	2^3	0 – open loop mode 1 – closed loop mode	0 8
4	2^4	0 – low pass filter off 1 – low pass filter on	0 16
5	2^5	0 – notch filter off 1 – notch filter on	0 32
6	2^6	reserviert	0
7	2^7	1 – real time signal processing (always on)	128
8	2^8	0 – simple current 1 – double current	0 256
9	2^9	0 – NanoX impossible 1 – NanoX possible	0 512
10	2^{10}	reserviert	0
11	2^{11}	actuator error (wrong actuator type)	2048
12	2^{12}	internal memory error	4096
13	2^{13}	I ² C-error	8192
14	2^{14}	underload*	16384
15	2^{15}	overload*	32768

table: status register

* Definition for overload and underload:

If it's impossible to reach the desired position in CL mode because of too much load or defect of the actuator (actual position is lower than desired), the overload condition is given. If the actuator not able to contract to lower position (actual position is higher than desired), the underload condition is given.

8.9 error messages

When user sends wrong inputs to the device, the controller responds with a message like “error,x”

The error numbers x means:

- 1 error not specified
- 2 unknown command
- 3 parameter missing
- 4 parameter range exceeded
- 5 parameter count exceeded
- 6 parameter is locked / read only

9 handling


The actuator is connected to the **NV100/D_NET** using the “PIEZO” plug. To control the amplifier by USB, connect the device USB connector to the PC USB port by the supplied USB cable and start a terminal application on PC (see chapter communication).

After switch on, the device runs a self test procedure, while the indicator LED is lightning red. This takes about 3 s. While this, the controller reads the system specific parameters from actuators ID-chip and set up the controller. Is there no actuator, the LED stay red.

After the self test, the actuator moves to its initial position. In case of open loop mode to the desired voltage, in case of closed loop mode it moves to the set position.

General to the OL and CL modes:

In OL mode, the actuator is direct controlled by a voltage between -20 V and +130 V. The reached stroke (pos. 1 to pos. 4) can be seen on actuators calibration protocol. The stroke is non-linear and shows hysteresis and also dependent by the actuators load. The stroke decreases by higher loads. To reach a defined position, actuators with integrated position sensors are used. These generates in CL mode a lower stroke, the remain stroke is used as a reserve to reach the end positions (pos. 2 and pos. 3) also with a (allowed) load.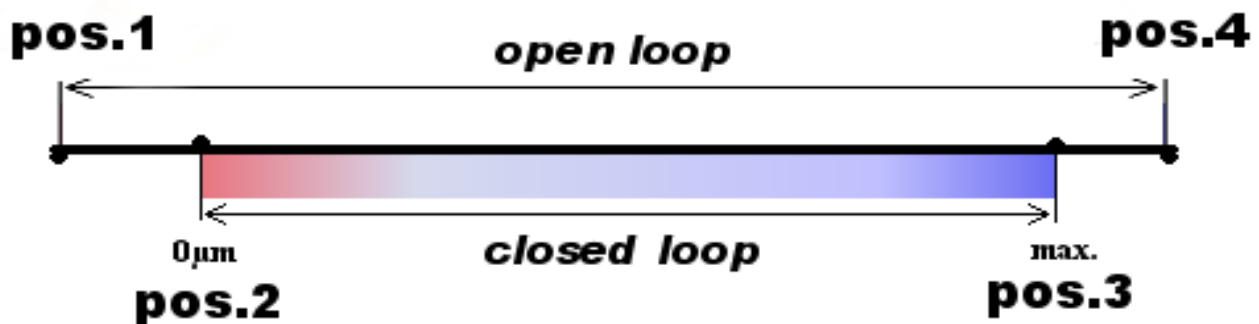

sketch 8: open loop / closed loop

When the desired position is not reached within 0.5 s, the device sets a flag in the status register (overload or underload flag). When the desired position is reached or a new position is set, the flags will be cleared and the overload or underload detection starts again. The state of the flags can be read with the status register.

10 controller adjustment

When any actuator made by **piezosystem jena** is connected to the **NV100/D_NET**, amplifier their specific values are read from the actuator's ID-chip. The Digital Signal Processor (DSP) of the amplifier is set with these values. These parameters were investigated in the **piezosystem jena** laboratory and ensure safe function of the actuator.

There are no k_p -, k_i -, and k_d -values for actuators without a measurement system because the closed loop mode is impossible in this configuration. Closed loop is automatically switched off.

To adapt the controller properties to your special application please start with the default values. At first please switch on the closed loop mode by the command **cl,1**. Change one parameter step by step and check the result. Maybe there is an additional external position sensor required.

In general, change the parameters in small steps and, depending on the actuator's reaction, slightly increase the step width. If the system begins to oscillate, switch off the closed loop (cl,0), then reset the last values entered!

Low dynamic applications need only an integral term in PID-controller, its set by parameter k_i , (k_p und k_d can be set 0). For filtering the set point, a slew rate limiter (command sr) is implemented. The limit is given in % / ms, it means, that a value of 1 %/ms would realize, that a jump off the full range (100 %) is converted to a ramp with 100 ms duration. The devices maximum slew rate depends on maximum output current and the actuators capacity. Independent from the sr value, this maximum slew rate cannot be exceeded.

To filter higher frequency parts on the input signal (in region to the resonant frequency or higher), a low pass filter 4th order is implemented. It can be activated or deactivated by the command lpon,1/0. The cut of frequency can be set by command lpf,value.

11 troubleshooting

Please check all cables and connections first if the system is not working properly.

error	possible solution
LED lights red	Check the connection between actuator and amplifier. Check the actuator cable for damage.
LED blinks red	NanoX is not possible on this device.
actuator oscillates in resonant frequency in closed loop	Check the PID-controller adjustments, reduce the ki-value, Reduce the kd-value.
no communication by USB is possible	<ul style="list-style-type: none"> - plug out and re plug in the USB cable, restart or reconnect the terminal application - check the device manager for a loaded USB driver - check the interface settings: - 115200 baud, 8 bits, 1 stop bit, no parity, software handshake (if applicable) Regardless the USB interface, the correct baud rate is necessary to set.
IP-Search did not find the device	<ul style="list-style-type: none"> - check the network connection cable - check the power supply (at powered device the led must light up) - check the firewall on PC and router (UDP 30718) - in the configuration „6 Security“ the setting „Port 77FE“ must set to “enabled” (use the browser to activate it)
telnet get the message: „the command ‚telnet‘ is wrong...“	<ul style="list-style-type: none"> - activate „Telnet Client“ under “Windows Features”
telnet don't get a connection, although the device is found in IP Search	<ul style="list-style-type: none"> - There is a connection with another program or from another computer to the device (it is only one connection possible) - The computer and the device are in different subnets <ul style="list-style-type: none"> o Set a static IP in the same subnet or o Change the subnet address on computer o Solve the problem with a direct connection to the PC („IP over DHCP“ must checked) - Check enabled firewalls (TCP 23) - In the configuration „1 Channel 1“ the „port“ is changed, check the setting

table: possible error causes

12 your notes